

# 3WAVES ELECTRA LITE XR-4 INSTALL

## Supplies Required

- Phillips screwdriver (provided)
- Cable guide and electrode holder (provided)



Ensure your Varjo XR-4 headset is powered off and unplugged before installation. Failure to accomplish this step may result in increased risk to the user or hardware.

1



- Remove the rear comfort pad from the XR-4.

2



- Align the 3WAVES Electra unit with three mounting holes in the rear of the headset. Use the provided screws to secure it into place.

3




- Reinstall the rear comfort pad in the the XR-4.

3



- Plug the forehead electrode lead (arrow) into the side of the Electra.

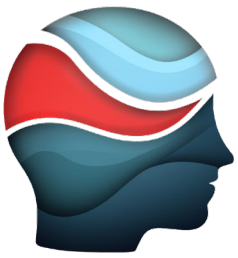

# 3WAVES ELECTRA XR-4 INSTALL GUIDE

## Supplies Required

- Phillips screwdriver (provided)
- Cable guide and wire receptacle (provided)



Ensure your Varjo XR-4 headset is powered off and unplugged before installation. Failure to accomplish this step may result in increased risk to the user or hardware.

5



- Remove the adhesive covers on the Cable Guide (A) and Wire Receptacle (B). Adhere them on headset as shown to secure front electrode lead.